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pH ELECTRODE FOR LOW CONDUCTIVITY WATER AND NON-AQUEOUS SOLVENTS INSTRUCTION MANUAL

In order to ensure correct usage of the pH electrode, be sure to read these operating instructions carefully before usage.

1. Specifications and the Electrode


This pH electrode has improved responsiveness in low conductivity water and non-aqueous solvents. It can be used with the F-10, M-10 and F-20 series of pH meters. The 6377 is a 3-in-1 electrode that contains a

glass electrode, reference electrode and temperature compensation electrode within a single body. The 1076A is a glass electrode. Refer to the diagrams below to verify the specifications for your electrode.

Type	6377-10D		1076A-10C	
Features	A 3-in-1 electrode with superior responsiveness in low conductivity, low bufferability aqueous solutions and non-aqueous solvents such as alcohol. It has a moveable sleeve type of liquid junction. The electrode can also be used for more general applications.		A glass electrode with superior responsiveness in low conductivity, low bufferability aqueous solutions and non-aqueous solvents such as alcohol. A separate reference electrode and temperature compensation electrode are required. The electrode can also be used for more general applications.	
Specifications	pH range Temperature range Internal electrode Internal solution for reference electrode Liquid junction Material of the parts that come in contact with the liquid pH meters supported	pH0~14 0~60°C Ag/AgCl 3.33 mol/L KCl solution (Cat. No. 300) Moveable sleeve Glass F-10, D-10, M-10, F-20 series	pH range Temperature range Internal electrode Reference electrodes supported Temp. electrodes supported Material of the parts that come in contact with the liquid pH meters supported	pH0~14 0~100°C Ag/AgCl 2060A, 2660A, 2565A 4163 Glass F-10, M-10, F-20 series
The electrode				

2. pH Meter Connection



- 1) Push the electrode connector as far as it will go onto the BNC terminal pin of the pH meter. Then rotate it to the right until it locks into place.
- 2) Cover the connector with the protective cover.
- 3) Push the temperature connector (for the 1076A, this is the connector for the corresponding temperature compensation electrode) as far as it will go onto the T terminal of the pH meter.
- 4) For the 1076A, attach the connector for the corresponding reference electrode to the R terminal of the pH meter.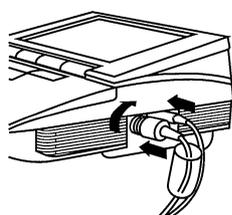

3. Procedures for First Time Usage or After Long-Term Storage


For the 1076A, perform these operations on the reference electrode.

- 1) Remove the rubber stopper from the internal solution filling opening and use a syringe to remove the internal solution.
- 2) Add new internal solution.
- 3) For the 6377 and sleeve type reference electrodes 2660A and 2565A, move the sleeve as shown in the diagram below. After checking the outflow of the internal solution, lightly pull the sleeve in the downward direction.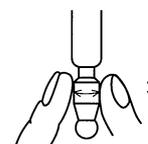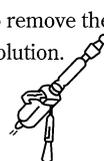


4. Measurement (Calibrations)


- 1) If there is not enough internal solution, remove the rubber stopper for the internal solution filling opening and add more internal solution.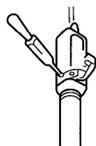

- 2) Wash the electrode more than 3 times with pure (deionized) water. Lightly wipe away any water droplets with tissue paper.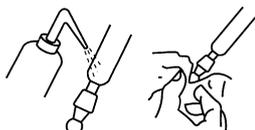

- 3) Immerse the electrode in the sample liquid. Be sure to keep the surface level of the internal solution of the electrode above the surface level of the sample liquid. Also be sure to remove the rubber stopper for the internal solution filling opening when taking measurements.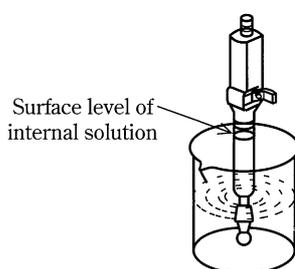


5. Precautions

<for both 6377 and 1076A>


- 1) Be sure to use Cat. No. 300 internal solution in the reference electrode. If you must use some other internal solution, use a 1076A glass electrode and 2565A double junction reference electrode combination. Place the appropriate internal solution in the outer chamber of the 2565A.
- 2) This electrode may be used to take measurements of alcoholic solutions which have very similar properties to water, and of non-aqueous solvents that contain a large proportion of water (greater than 5%). However, the responsiveness and stability is not as good as for aqueous solutions.
- 3) The electrode connectors require a high degree of insulation. Do not let them get wet or touch them with your hands.
- 4) When a high degree of accuracy is required, after calibration, wash the detection apparatus more than 3 times with pure (deionized) water and soak it for 10 minutes more in pure (deionized) water before taking measurements

<6377>

- 1) If there are any bubbles on the inside of the liquid junction opening of the sleeve liquid junction, or in the thin part of the glass pipe, shake the electrode lightly up and down to free the bubbles.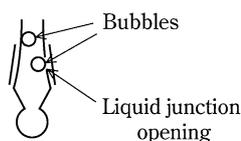

- 2) If the sleeve liquid junction becomes hardened with precipitated crystals from the internal solution and will not move, soak it in warm water of about 30 to 40 degrees to dissolve the precipitate. Move the sleeve liquid junction taking care not to damage it. Check the outflow of the internal solution before usage. (Submitting the sleeve liquid junction to an ultrasonic cleaner is also effective.)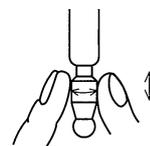

- 3) Press the sleeve lightly down. If you push it too strongly, it will cause a large increase in the liquid junction potential and give rise to measurement errors. If the sleeve is not closed tightly enough then the internal solution may seep out too rapidly, or the sample solution will seep in. If the sample solution seeps in, it will cause a large measurement error, and the internal solution will have to be replaced.

6. Regular Maintenance

- 1) If the electrode is used for long periods, the internal solution in the reference electrode may become contaminated by the sample liquid, or become diluted. Perform the operations described in step 3. Procedures for First Time Usage or After Long-Term Storage once every 3 to 6 months.
- 2) If the inside of the protective cap becomes dry, wash it with pure (deionized) water. After removing any excess water, supplement it with enough pure (deionized) water to soak the sponge.
- 3) If the response time of the electrode or its reproducibility has fallen, wash the electrode using the procedures described below. Rinse it well with pure (deionized) water and soak only the glass membrane for several hours in pure (deionized) water before taking measurements.

Ordinary dirt	: Wipe with gauze soaked in a neutral detergent.
Oily dirt	: Wipe with gauze soaked in an appropriate organic solvent such as alcohol.
Dirt containing inorganic components	: Soak for approximately 2 hours in 1 mol/L hydrochloric acid. Be sure not to immerse the liquid junction in the hydrochloric acid.